



Kendall™ Disposable Tissue Grinder Systems

Single Use System

- No cleanup or sterilization required
- Grinder is completely disposable after use

Protective Sleeve

- Contains any aerosolization created during grinding

Hollow Pestle

- Ice can be added to pestle core to keep sample cool

Ergonomic Handle

- Easy grinding

Built in Transportation System

- No transfer of specimen required for transportation or storage



Disposable Tissue Grinder Systems



Item Number	Description	Case Quantity
3505SA	Small Tissue Grinder with 15mL tube	10
3500SA	Large Tissue Grinder with 50mL tube	10

Instruction For Use



Place tissue in tube on side wall. Fluid can be added if desired.



Hold tube sideways and insert pestle, making sure tissue is pushed to bottom of tube.



Hold tube through sheath and grind back and forth through 90's (degrees).



To rinse pestle tip, back pestle out of tube 1/4 inch (6.3mm) and shake side to side.



After grinding discard pestle in appropriate manner.



Cap tube and store or transport.



COVIDIEN, COVIDIEN with logo, Covidien logo and *positive results for life* are U.S. and/or internationally registered trademarks of Covidien AG. TM trademarks of a Covidien company.
©2009 Covidien.
H6498 5M 1109

For Ordering Information Contact CHARTECH INC.
1-877-658-1849 Fax 770-442-2843
sales@chartech.net